

PICO Protein Detection Trial Kit USER MANUAL



Actome GmbH Georges-Köhler-Allee 103 79110 Freiburg Germany

01 Kit Contents and Storage	3
02 Intended Use	3
03 Safety Information	4
04 Quality Control	4
05 Introduction	5
06 Scientific Background of a CLC Experiment 06.1 Explanation of the Bell-shaped Curve 06.2 Using CLC Experiments for Relative Quantification in Different Biological Samples	6 6 7
07 Preparation for PICO Assay	9
07.1 General Remarks 07.2 Plate Layout for Setup of Binding Reaction and dPCR Pre-Dilution	9 10
08 Equipment and Reagents to be Supplied by User	11
09 Protocol of Protein Detection Trial Kit	12
09.1 Buffer Preparation	12
09.2 Preparation of the Dilution Series	13
09.3 Binding Reaction	13
09.4 Pre-dilution and Digital PCR	15
10 Evaluation	18
10.1 Raw Couplex Calculation using AMULATOR	18
10.2 Couplex Count Processing for Relative Quantification (RQ)	19
10.2.1 Lambda Normalization	19
10.2.2 ABC Correction	21
10.3 Relative Quantification (RQ)	22
11 Troubleshooting Guide	25
12 Ordering Information	26



01 Kit Contents and Storage

PICO Protein Detection Trial Kit				
Catalog Number:		PICO-000090		
Pouch - store at 4°C	pcs	Item code		
Additive T	1	А		
Additive C	1	В		
Bovine serum albumin	1	С		
Prelabeled Trastuzumab antibody $(4.82 \times 10^9 \text{ cp/}\mu\text{l})$	1	D		
Prelabeled Pertuzumab antibody $(4.82 \times 10^9 \text{ cp}/\mu \text{l})$	1	Е		
Pouch - store at -20°C				
Additive L	1	F		
Coupling dPCR Mix	1	G		
Recombinant human ErbB2/HER2 protein (0.529 µmol/l)	1	Н		
PICO BL Probe	1	Ι		
PICO P8 Probe	1	J		

The PICO Protein Detection Trial Kit is shipped at room temperature. Upon arrival, it is important to ensure that the products are immediately transferred to optimal storage conditions. Please note the different storage temperatures for the components of the PICO Protein Detection Trial Kit. Date of expiry is stated on the label of each package.

02 Intended Use

The PICO Protein Detection Trial Kit is intended for molecular biology applications only. The product is not intended for the diagnosis, prevention, or treatment of a disease. All due care and attention should be exercised in the handling of the product. We recommend the users of the Actome products to adhere to the national safety guidelines.



03 Safety Information

When working with chemicals, always wear a suitable lab coat, disposable gloves, and protective goggles. For more information, please consult the appropriate material safety data sheets (MSDS). These are available at: www.actome.de/resources/downloads

In case of reagent spillage, absorb the spilled material, dispose of it accordingly and clean with suitable laboratory detergent and water. If the spilled liquid contains potentially infectious agents, clean the affected area first with laboratory detergent and water, and then with 1% (w/v) sodium hypochlorite.

04 Quality Control

Each lot of PICO Protein Detection Trial Kit is tested against predetermined specifications to ensure consistent product quality.



05 Introduction

Protein Interaction Coupling (PICO) is a highly sensitive immunoassay for the detection of proteins, protein interactions, and post-translational modifications using QIAGEN's QIAcuity Digital PCR System. The **PICO Protein Detection Trial Kit** is an ideal tool for researchers who want to become familiar with the PICO assay. It includes all the necessary reagents to perform one PICO measurement, using the recombinant human ErbB2/HER2 protein as the target, along with two pre-labeled antibodies (trastuzumab and pertuzumab) that are directed against the target. The Trial Kit highlights the unique features of PICO, including ultra-high sensitivity, zero background, wash-free workflow, and high specificity. With the PICO Protein Detection Trial Kit, you'll gain a deep understanding of the core PICO technology, experimental set-up, and data analysis using the AMULATOR. Additionally, the kit comes with a detailed protocol for setting up a calibration curve experiment (CLC), which involves using a fixed antibody concentration and a dilution series of the sample input to demonstrate the relative quantification (RQ) evaluation of the assay (**Figure 1**).

If you would like to design your own PICO assay consider the following instructions:

The PICO assays can be performed using two, three, or four different labeled antibodies, depending on how many targets you would like to measure. A target can be a protein, a protein interaction, or a post-translational modification. For the detection of an individual target two different labeled antibodies are required. In the case of a single protein detection the two antibodies must bind to non-overlapping epitopes. For the detection of a protein interaction one antibody per interacting partner is required. To detect a post-translational modification (PTM) a PTM-specific antibody and an antibody directed against the PTM-bearing protein is required. The molecular complex, consisting of the target and two bound antibodies (called 'couplex'), is the molecular detection unit of the PICO assay.

For the PICO assay the antibodies of your choice are labeled with **PICOglue Labels** (available in four colors: BL, P8, N6, or O7) using the **PICOglue Antibody Labeling** (gAL) Kit. Next, the **PICO Amplification Core Kit** is used to perform the PICO reactions. If the antibody solution contains any chemicals, such as BSA, azide, TRIS, or free primary amines, we recommend using the **PICO Antibody Purification Kit** prior to using the PICO gAL Kit.

Day	Steps	Timing	Hands-on	Stopping points & Storage
DAY 1	Step 1 – Preparation of PICO Assay	55 min	15 min	
	A. Preparation of Dilution Series of Protein	30 min	10 min	
	B. Setup of Antibody Mix	25 min	5 min	
	Step 2 – Calibration Curve	4 h	40 min	
	A. Binding Reaction	30 min + ON	10 min	overnight incubation (ON) at 4°C
DAY 2	B. Dilution of the Samples before dPCR	20 min	20 min	
	C. QIAcuity dPCR Run	3 h 10 min	10 min 🛛 🧧	overnight run possible
	Step 3 – Evaluation	35 min	25 min	
	A. Raw Couplex Calculation using AMULATOR	15 min	5 min	
	B. Relative Quantification	20 min	20 min	

Figure 1. The PICO Protein Detection Trial Kit workflow.

06 Scientific Background of a CLC Experiment

06.1 Explanation of the Bell-shaped Curve

The Protein Detection Trial Kit protocol includes a calibration curve (CLC) experiment to determine the optimal linear range between antibody and target concentration. In the protocol a dilution series of the biological sample with a constant antibody concentration is prepared. If the antigen concentration in the sample is high enough, the couplex counts derived from a CLC experiment can form a bell-shaped curve when plotted against the dilution series of the sample input (**Figure 2**).



Figure 2. Typical result of measuring a dilution series with PICO: a bell-shaped curve Boxplots representing the absolute number of couplexes detected at different dilution steps (from highest 'S#5' to lowest 'S#1' dilution). The pattern of the boxplots resembles a bell-shaped curve. Ratios between antigen and labeled antibodies are depicted at three different dilution steps (S#5, S#3, and S#1). Couplexes are highlighted with a yellow background. The central solid line within each box is the mean, and the bottom and top of each box represent the 25th and 75th percentiles, respectively. The whiskers represent the 10th and 90th percentiles. ABC - antibody control

Diluting the biological sample leads to a decreasing amount of antigen in the binding reaction (the mixture of biological sample and labeled antibodies). In **Figure 2** the absolute number of couplexes, measured from a 3-fold dilution series of recombinant HER2 protein using 4×10^{-11} M antibody concentration in the binding reaction, is depicted. The S#1 sample contains the highest antigen amount (1×10^{-9} M), while the S#5 sample has the lowest antigen amount (1.23×10^{-11} M). The <u>Antibody Control (ABC)</u> is a control sample containing only the labeled antibodies.

On the left side of the curve (low antigen side), the ratio of antibody to antigen is high, leading to decreased probability of couplex detection (S#5 to S#4 in **Figure 2**). The high dilution of the sample leads to a loss of couplex count in the dPCR, since the couplexes are diluted out and the compartments are occupied mainly by unbound labeled antibodies.

Increasing the antigen concentration leads to an increasing probability of couplex formation. In a certain range, this trend is linear until an equilibrium between antigen and antibodies is reached (approximately at S#3 in **Figure 2**). At this equilibrium, the highest amount of couplexes are measured.



On the right side of the curve (high antigen side), the further increase in antigen concentration leads to a decrease in couplex count (S#2 to S#1 in **Figure 2**). The increase of antigen amount decreases the chance of couplex formation, since the chance that the two different labeled antibodies bind to the same target is smaller. In a certain range this trend is linear until the number of partitions of the dPCR is too low for couplex detection.

The linear correlation between the couplex counts and the antigen amount at the low antigen side of the curve (S#5 to S#4 in **Figure 2**) can be used for relative quantification (RQ), to relatively compare the amount of target between different sample types. For this, the raw couplex counts are processed and the molar ratio between the couplex counts of two different samples is calculated (see 10.2 and 10.3). The normalization of sample input allows the comparison of data derived from different amounts of sample. However, for relative comparison of two samples, the same antibody concentration in the binding reaction has to be used. Additionally, for better comparability between experiments, we recommend to use the same batch of labeled antibodies.

06.2 Using CLC Experiments for Relative Quantification in Different Biological Samples

Careful planning of an CLC experiment is necessary to generate a dataset that can be used for relative quantification. See the different outcome scenarios for a CLC experiment in **Figure 3**.



Figure 3. Hypothetical outcome scenarios from a calibration curve (CLC) experiment.

Boxplots representing the absolute number of couplexes detected at different dilution steps (from highest 'S#5' to lowest 'S#1' dilution). (A) PICO data reflecting the low antigen side of the bell-shaped curve, indicated by increasing couplex count with increasing antigen concentration. (B) PICO data reflecting the high antigen side of the bell-shaped curve, indicated by decreasing couplex count with increasing antigen concentration. (C) PICO data with low couplex counts due to over diluted samples or low expression levels of the measured target. The central solid line within each box is the mean, and the bottom and top of each box represent the 25th and 75th percentiles, respectively. The whiskers represent the 10th and 90th percentiles. ABC - antibody control.

In general, biological samples with a high expected concentration of the measured target require less sample input in the binding reaction, while biological samples with a low concentration of the measured target require more sample input. For initial experiments we recommend starting from a cell lysate equivalent of 20,000 cells in the binding reaction and performing a 2-fold (low expressing target) or 3-fold (high-expressing target) dilution series. Since biological samples vary widely, these settings may need to be adjusted for each sample. In case of using recombinant protein material, we recommend to perform a dilution series where the concentration of the dilution step S#3 is in the same range as the antibody concentration used in the binding reaction (in most cases 4×10^{-11} M).

In an optimal case, a CLC experiment results in a curve as depicted in **Figure 3A**. In this example the antibody concentration in the binding reaction is 4×10^{-11} M. The S#1 sample



contains the highest protein amount $(4 \times 10^{-11} \text{ M})$ while S#5 has the lowest protein amount $(4.94 \times 10^{-11} \text{ M})$. Increasing the sample input leads to an increase in couplex count, indicating that the measured data is located on the low antigen side of the bell-shaped curve. This CLC pattern is observed for most experimental setups, especially if the expression of the target molecule is rather low. Data points derived from dilutions in the linear range of this curve can be used for RQ. We recommend not to use the data from the highest measured couplex count to ensure that only data from the linear range is used.

In another scenario, a CLC experiment might result in a curve as depicted in **Figure 3B**. In this example the antibody concentration in the binding reaction is 4×10^{-11} M. The S#1 sample contains the highest protein amount (3×10^{-8} M) while S#5 has the lowest protein amount (1.23×10^{-10} M). Increasing the sample input leads to a decrease in couplex count, indicating that the measured data is located on the high antigen side of the bell-shaped curve. This CLC pattern is observed for experimental setups with a high expression level of the target. In this case, we recommend repeating a CLC experiment with higher dilution factors for the biological samples so that the results reflect couplex counts as seen in **Figure 3A**. However, this scenario will be observed only with very highly expressed targets.

In a third scenario, a CLC experiment might result in a curve as depicted in **Figure 3C**. In this example the antibody concentration in the binding reaction is 4×10^{-11} M. The S#1 sample contains the highest protein amount $(1.27 \times 10^{-12} \text{ M})$ while S#5 has the lowest protein amount $(1.57 \times 10^{-14} \text{ M})$. Very low couplex counts are measured for all dilution steps. In this case, the expression level of the target might be too low for the chosen experimental setup. We recommend increasing the overall sample input (e.g. using a higher amount of lysed cells) or decreasing the dilution factors for the biological sample.

The CLC experiment described in the PICO Protein Detection Trial Kit protocol results in a curve as shown in **Figure 2**. It demonstrates the full bell-shaped curve at the different antigen-to-antibodies ratios.



07 Preparation for PICO Assay

07.1 General Remarks

- The PICO assay is sensitive to variations in pipetting. Make sure all pipettes are calibrated for reliable liquid handling. The dilution factor for dPCR is usually very high and it is important to use highest pipetting precision available to ensure reliable quantitative results (see recommended equipment and devices in 08 Equipment and Reagents to be Supplied by User).
- Use gentle mixing by pipetting up and down. This helps to reduce spill-overs and loss of material.
- Change gloves regularly during laboratory work to avoid contamination. Contamination can be detected if the NTC sample shows positive partitions in any color.
- To ensure the highest molecular sensitivity and reproducibility of PICO we recommend using four technical replicates of each sample.
- In addition to the samples and the <u>non template control (NTC for dPCR)</u>, <u>antibody control</u> (ABC) reactions are also set up. An ABC reaction contains the corresponding antibody mix without the sample. In these samples, the two different labeled antibodies colocalize in the compartments only according to Poisson statistics. Therefore, the couplex count in the ABC should be zero. However, due to offsets in the dPCR (e.g. signal dropouts or incorrect clustering) the ABC value may deviate from zero. Thus, the AMULATOR performs an automatic ABC normalization by default.
- The average number of targets per partition is called lambda. Lambda is also a statistical parameter for the underlying Poisson distribution. A higher lambda value means higher assay sensitivity, but also a larger standard deviation based on Poisson statistics and a loss of signal due to dropouts, since several templates have to be amplified in one partition. For the PICO assay, we recommend a lambda of 0.15, however any lambda values between 0.01 and 0.6 can be used for PICO.
- To reach an optimal lambda for the PICO assay, the samples must be diluted prior to the dPCR. The dilution of the samples prior to dPCR disrupts the equilibrium binding conditions and induces dissociation of the antibodies from their target. After the samples have been diluted, make sure to work without interruption until the QIAcuity Nanoplate 26k 24-well is loaded into the QIAcuity.





07.2 Plate Layout for Setup of Binding Reaction and dPCR Pre-Dilution



The binding reaction setup and the dPCR pre-dilution for a typical PICO assay can be seen in **Figure 4**. First the dilution series of the sample is combined with the antibody mix and incubated overnight (binding reaction). We recommend preparing three technical replicates for the ABC and four technical replicates for each sample from the binding assay. The following day the binding reaction is diluted for the dPCR reaction. First, a 96-well plate is prepared by adding the calculated amount of PBS to columns 1 -3 and 5 - 7. In column 10 - 12 41 µl of Master Mix is added. Then, the binding reaction is diluted 1:10 using PBS and mixed by pipetting 30 times ('pre-dilution of the binding reaction'). From each diluted binding reaction sample, 1 µl is transferred into the corresponding wells of the first dilution step (DS1) and the dilution is mixed by pipetting 30 times. Then, 1 µl of DS1 is transferred to the corresponding wells of the second dilution step (DS2) and the dilution is mixed by pipetting 30 times. Finally 1 µl of DS2 is transferred to the wells containing the Master Mix. After mixing by pipetting three times, 40 µl of the Master Mix containing the diluted sample are transferred to a QIAcuity Nanoplate 26k 24-well and dPCR is performed. Perform the dilution steps as fast as possible since the dilution disrupts the equilibrium binding conditions and induces antibody dissociation.

08 Equipment and Reagents to be Supplied by User

Devices/Equipment

- QIAcuity Digital PCR System (QIAGEN, Cat.#: 911001)
- Table-top mini centrifuge for quick spins (~1,000 rcf)
- Plate centrifuge (e.g. Megafuge 8, Thermo Fisher, Cat.#: 75007210)
- Vortex mixer
- Ultrasonic bath
- Multichannel pipette, 8-channel $(1 10 \mu l)$
- Multichannel pipette, 8-channel $(10 100 \mu l)$
- Regular 1-channel pipettes (1 1,000 µl)
- Electronic Multichannel pipette (INTEGRA, Cat.#: 4722), 12-channel, 5 – 12.5 μl (optional)

Consumables

- QIAcuity Nanoplate 26k 24-well (QIAGEN, Cat.#: 250001)
- 2 PCR microplates, 96 well, Polypropylen, V-bottom or similar
- Sealing foil Adhesive Film (e.g. Thermo Fisher, Cat.#: 10696771)
- 12.5 µl GRIPTIP, Sterile, Filter 5 Racks of 384 Tips, long; for electronic multichannel pipette (INTEGRA, Cat.#: 3405) (optional)
- 1.5 ml reaction tubes
- 0.5 ml reaction tubes
- 10 µl, 200 µl, 1,000 µl standard pipette tips

Chemicals and Kits

- QIAcuity Probe PCR Kit (QIAGEN, Cat.#: 250101)
- Phosphate-Buffered Solution (PBS), without calcium or magnesium ions (e.g. Thermo Fisher Scientific, Cat.#: 12037539)
- cOmplete Protease[™] Inhibitor Cocktail (Roche, Cat.#: 04693132001)
- Ultrapure water (e.g. Thermo Fisher Scientific, Cat.#: 15667708)



Notes

09 Protocol of Protein Detection Trial Kit

Note: The protocol is valid for performing a single PICO assay using the PICO Protein Detection Trial Kit.

09.1 Buffer Preparation

1. Prepare the chemicals and buffers as listed below. Prepare the buffers directly before they are used.

Additive C (5x stock) Add 500 μl PBS Easily soluble, stable for 1 week at 4°C.

BSA (5x stock) Add 400 µl PBS Stable for 3 days at 4°C

EDTA-free Protease Inhibitor Cocktail (PIC), (25x stock) 1 tablet of cOmplete Protease[™] Inhibitor Cocktail 2 ml PBS Stable for 12 weeks at -20°C

Lysis Buffer Stock (LB-Stock), (2x stock) 200 µl Additive T 400 µl Additive C 80 µl PIC 200 µl Additive L 120 µl PBS Stable for 3 days at 4°C

Lysis Buffer (LB) 300 µl LB-Stock 300 µl PBS Stable for 3 days at 4°C

Control Buffer (CB) 250 µl LB-Stock 100 µl BSA 150 µl PBS Prepare the Control Buffer fresh



09.2 Preparation of the Dilution Series

- 2. Thaw a vial of the supplied recombinant human ErbB2/HER2 protein and spin it down (~1,000 rcf, 5 s).
- 3. Prepare the dilution series of the protein in LB:

Start with a 100-fold dilution (sample S#1). Next, dilute sample S#1 60-fold (total dilution factor: 6,000) to generate sample S#2. Dilute sample S#2 2-fold (total dilution factor: 12,000) to generate sample S#3. Dilute sample S#3 2-fold (total dilution factor: 24,000) to generate sample S#4. Finally, dilute sample S#4 10-fold (total dilution factor: 240,000) to generate sample S#5. Prepare a volume of at least 30 µl for each sample. If required, plan the dilution series using the <u>PICO</u> <u>Calculator</u> (step 1 in the spreadsheet).

09.3 Binding Reaction

- 4. Prepare the antibody mix (ABX) in a total volume of 100 μ l. Calculate the volume of antibody stock, LB-Stock, and PBS added to the ABX by following steps 5 to 7. Alternatively, use the <u>PICO Calculator</u> (step 2 in the spreadsheet) for the calculations and proceed directly to step 8. Use an antibody concentration of 4 × 10⁻¹¹ M and a dilution factor of 2 in the binding reaction.
- 5. To calculate the volume of each antibody stock added to the ABX, the concentration of the antibody stocks must be converted from copies per μ l to molar concentration.

$$C_{M} = C_{C} \times 10^{6} / N_{A}$$

 $C_{_M}$: Conc. of antibody stock [M]

 C_c : Conc. of antibody stock $[cp/\mu l]$

This concentration is determined via dPCR in the ABQC (see user manual aCALL Kit)

 N_{A} : Avogadro constant (6.02214076 × $10^{23} mol^{-1}$)

Example: The concentration of the supplied antibody stock solutions (C_c) is 4.82×10^9 cp/µl. The calculated molar concentration is (8×10^{-9} M).

6. Calculate the volume of each antibody $(V_{AB1} \text{ or } V_{AB2})$ used in the ABX based on the following calculation.

If V_{ABI-2} are too low for precise pipetting, prepare a pre-dilution of the antibody stock using LB. In this case, the concentration of the pre-dilution of the antibody (DF_{AB}) has to be considered in the calculation.

$$V_{_{AB}} = V_{_{ABX}} \times C_{_{BRM}} \times DF_{_{BR}} \times DF_{_{AB}}/C_{_{M}}$$

 V_{AB} : Volume of antibody stock (or dilution) added to the ABX V_{ABX} : Total volume of ABX; we recommend at least 60 µl C_{BRM} : Conc. of antibody in the binding reaction [M];

we recommend to use 4 \times 10⁻¹¹ M

 DF_{BR} : Dilution factor introduced by pipetting of binding reaction DF_{AB} : Dilution factor introduced by pre-dilution of antibody stock ; use '1' if no predilution is performed

Example: The molar concentration of the antibody stock is 4.82×10^{-9} M. We choose 100 µl as the total volume of the ABX and an antibody concentration of 4×10^{-11} M in the binding reaction. The dilution factor introduced by the binding reaction setup is 2 in our case (2 µl of the ABX is combined with 2 µl of the cell lysate). The calculated volume of each antibody stock added to the ABX is 1 µl.

7. Prepare the ABX using LB-Stock. Add the calculated volume of the two antibody stock solutions or prepared dilutions and fill up the remaining volume with PBS.

If a high volume of antibodies has to be used, the amount of PBS added to the ABX might be negative, in this case use LB to fill up the whole ABX volume.

Example: The total volume of the ABX is 100 μ l. Add 50 μ l of LB-Stock in a low protein binding tube. The volume of each antibody added to the ABX is 1 μ l. Add 48 μ l PBS to fill up the volume to 100 μ l.

- Set up the binding reaction in a 96-well PCR microplate (v-bottom) by mixing 2 μl of ABX with 2 μl of S#1-S#5 (Figure 4). Make four technical replicates for each dilution step.
 Due to the small volume of the binding reaction, do not mix the ABX with the biological sample by pipetting.
- 9. For the ABC, combine 2 μ l of CB with 2 μ l of ABX and make three technical replicates.
- 10. For the NTC add 4 μ l of LB to one dedicated well.
- 11. Seal the plate (without the lid) with an adhesive foil. Sonicate the plate at full power for 1 min in an ultrasonic bath. Centrifuge the plate to collect the liquid at the bottom (~1,000 rcf, 30 s) and incubate at 4°C overnight. The incubation time can be varied between 12-24 hours.

09.4 Pre-dilution and Digital PCR

12. Prepare the Master Mix for 24 samples (see amounts below). Vortex for 10 s and spin down (~1,000 rcf, 5 s). Store the Master Mix at 4°C.

Master Mix:

697 μl Ultrapure water 284 μl QIAcuity Probe Master Mix 45 μl PICO P8 Probe 45 μl PICO BL Probe 36 μl Coupling dPCR Mix

- 13. To reach the optimal lambda range for dPCR, the samples need to be diluted first. We recommend aiming for a lambda of 0.15 and performing a 10x pre-dilution of the binding reaction due to the low volume. The dilution factor introduced by the Master Mix is 42. To calculate the necessary dilution factor for a lambda of 0.15, follow steps 14 to 17 or use the <u>PICO Calculator</u> (step 3 in the spreadsheet) for the calculations and proceed directly to step 18.
- 14. Convert the antibody concentration, used in the binding reaction, from molar concentration to copies per μl.

$$\begin{split} C_{BRC} &= C_{BRM} \times N_A / 10^6 \\ C_{BRC} : Conc. of antibody in the binding reaction [cp/µl] \\ C_{BRM} : Conc. of antibody in the binding reaction [M] \\ N_A : Avogadro constant (6.02214076 \times 10^{23} mol^{-1}) \end{split}$$

Example: The concentration of antibodies in the binding reaction is 4×10^{-11} M. The concentration in cp/µl is 2.4×10^{7} .

15. Calculate the targeted antibody concentration for the dPCR.

 $C_d = \lambda / V_p$

 C_d : Targeted conc. of antibodies for the dPCR [cp/µl]

 λ : Average number of targets per partition: we recommend 0.15

 V_n : Volume of a partition [µl]: 0.00078 µl

Example: We choose a lambda of 0.15. The volume of one partition of a QIAcuity Nanoplate 26k 24-well is 0.00078 μ l. Thus, the targeted concentration for the dPCR is 192.31 cp/ μ l.

16. Calculate the dilution factor necessary to reach an optimal lambda range for the dPCR (DF_d). Based on the small volume of the binding reaction, we recommend to perform a pre-dilution of the binding reaction. This, together with the dilution introduced by the volume of the Master Mix has to be considered in the calculation of the DF_d.

$$DF_d = C_{BRC} / (C_d \times DF_{pre} \times DF_{MM})$$

 DF_d : Dilution factor necessary to reach an optimal lambda DF_{pre} : Dilution factor of pre-dilution; we recommend 10 DF_{MM} : Dilution factor introduced by Master Mix; usually 42

Example: The concentration of antibodies in the binding reaction is 2.4×10^7 copies/µl and the targeted concentration for the dPCR is 192.31 copies/µl. Considering a 10x pre-dilution of the binding reaction and the dilution introduced by the volume of the Master Mix, the dilution factor is 298.24 (which can be rounded to 298). This means 1 µl of the 10x pre-diluted sample is diluted in 297 µl PBS.

17. Normally, the necessary dilution volume to reach the optimal lambda range is very high. For easier handling we recommend to perform a serial dilution using the following calculation.

Please consider that the V_d is rounded for easier handling of the pipetting.

$$V_d = (((DF_d)^{\frac{1}{\#DS}}) \times V - V_c)$$

 V_d : Volume [µl] of PBS required for each dilution step #DS: Number of dilution steps V_c : Carry-over volume of previous dilution; usually 1 µl

Example: The dilution factor corrected for the 10x pre-dilution and the dilution introduced by the volume of the Master Mix is 298.24 (which can be rounded to 298). To prevent using a large dilution volume we chose a serial dilution including two dilution steps (#DS = 2) and a carry-over volume of the previous dilution of 1 µl. This results in a volume of 16.27 µl PBS which is rounded to 16 µl. This means, 1 µl of the 10x pre-diluted binding reaction samples are transferred to 16 µl of PBS for the first dilution step and 1 µl of this dilution step will be transferred to 16 µl of PBS for the second dilution step.

18. Prepare a 96-well plate with columns 1 - 3 and 5 - 7 containing the calculated amount of PBS for the dilution steps and columns 10 - 12 containing 41 μ l of Master Mix (**Figure 4**).

- 19. Remove the adhesive foil from the plate containing the binding reaction carefully. Add 36 µl of PBS to the wells (represents the first 10x pre-dilution). Mix thoroughly by pipetting up and down 30 times.
 Important: Perform the following steps without unnecessary breaks as the dilution disrupts the equilibrium binding conditions and induces antibody dissociation.
- 20. Transfer the predetermined carry-over volume from each diluted sample into the corresponding wells (columns 1 3) of the dilution plate. Mix by pipetting up and down 30 times (Figure 4). This dilution step is defined as DS1.
- 21. Transfer the predetermined carry-over volume from the first dilution step (DS1) into the next set of corresponding wells (columns 5 7). Mix by pipetting up and down 30 times (Figure 4). This dilution step is defined as DS2.
- 22. Transfer 1 μ l from the last dilution step (DS2) into the corresponding wells containing 41 μ l Master Mix (columns 10 12). Mix by pipetting up and down 30 times (**Figure 4**).
- 23. Finally transfer 40 μl of the Master Mix containing the diluted samples into a QIAcuity Nanoplate 26k 24-well. Seal the plate according to the QIAcuity user manual protocol and insert the plate into the QIAcuity dPCR System. Run the dPCR program on the QIAcuity according to the instructions of the manufacturer using the following dPCR settings:

Priming

QIAGEN Standard Priming Profile

PCR conditions

Hot-start	95°C for 2 min
Cycling 40 times	
Denaturing	95°C for 15 s
Annealing	58°C for 30 s

Imaging conditions

PICO P8 Probe:FAM green channel, 500 ms integration time, gain 6PICO BL Probe:HEX yellow channel , 400 ms integration time, gain 6



10 Evaluation

10.1 Raw Couplex Calculation using AMULATOR

24. After the dPCR run, download the raw dPCR data (RFU values) from the QIAcuity Software Suite. Select your plate in the 'Plates Overview', click 'Analyze' and select all wells. In the 'Select targets' drop down window click 'Select All' and press 'Show results'. Click 'Export to CSV...' and select 'RFU values'. The download generates a zip file containing csv files (one for each fluorescent color). Unzip the exported CSV file.

Please control the raw data and images of the plate in the QIAcuity Software Suite to ensure that the data is valid and the experimental setup was performed correctly.

25. Prepare the <u>Sample Definition File</u>. The spreadsheet contains two sheets: 'Layout' and 'Settings'. In the 'Layout' sheet fill in the name of the samples ('SampleName') to the corresponding wells. Replicates are defined by having the same sample name. In the 'ABC' column mark the ABC samples with 'true'. In the 'ExperimentalGroup' column, group together the samples with their corresponding ABC control by using the same number. If the same ABX is used for all samples, use the same number. The NTC sample is in its own group.

Important: Sample names should contain letters (e.g. sample #1). In case of invalid wells (wells containing no signals) the corresponding wells in the sample definition file as well as the according table line have to be deleted to create a continuous table.

- 26. In the 'Settings' sheet of the sample definition file, assign the labeled antibodies to the corresponding fluorescent color according to the PICO labels used (FAM-P8, HEX-BL).
- 27. Create a zip file containing the RFU files and the sample definition file. The data is now ready to be uploaded to the AMULATOR software.
- 28. Access to the <u>AMULATOR</u> software is provided on the website. The AMULATOR user manual is available <u>here</u>.
- 29. First-time users require an initial registration to our server. Registration provides a personal cloud-based storage space, where you can store and (re)analyze your uploaded PICO datasets (up to 15 files). The files are accessible only by you.
- 30. Upload the zip file using the ten-character token provided with the kit. The uploaded files are organized in a table. By clicking on the delete sign you can permanently delete them (after confirmation).
- 31. Mark the 'Add' checkbox next to the dataset you want to analyze. Click on the filename to initiate the data analysis. A red spinning box appears indicating that the data is being processed. When the analysis is finished, the box turns green.

Important: Wait until the analysis is finished (box turns green) before continuing.

- 32. The AMULATOR creates two graphs. The boxplots in the couplex chart display the number of couplexes for each sample, including the ABC controls. The boxplots in the lambda charts represent the lambda values of each antibody in each well. In the couplex chart each individual boxplot can be arranged by drag and drop. To delete a single datapoint or an entire boxplot hover over the data and double click on the popup text. To apply ABC subtraction, use uniform Y axis, or to display the standard statistics of normality use the checkboxes below the charts.
- 33. To download the results click on the 'Download PICO results' button. To download the graphs click on the 'Download couplex_charts' or 'Download lambda_charts' button.
- 10.2 Couplex Count Processing for Relative Quantification (RQ)

Note: The experimental setup of the PICO Protein Detection Trial Kit allow us to perform RQ and relatively compare the amount of recombinant human ErbB2/HER2 protein in the different sample dilutions on the low antigen side (see *02 Scientific Background of a CLC Experiment*). To demonstrate the bell-shaped curve behavior of the PICO assay, the dilution series in the PICO Protein Detection Trial Kit is set in a manner that S#2 to S#5 are on the low antigen side, while S#1 is on the high antigen side. Therefore, we can use S#2 (6000-fold diluted sample), S#3 (12,000-fold diluted sample), S#4 (24,000-fold diluted sample) and S#5 (240,000-fold diluted sample) for RQ. Sample S#5 is over-diluted to demonstrate the lower detection limit of the assay and it is a statistically non-significant signal above zero. This will be demonstrated mathematically at *section 10.3 step 45*.

- 34. The raw couplexes from the AMULATOR are used as an input for RQ. Download the raw couplex count <u>without ABC correction</u> from the AMULATOR software (see *10.1 step 32*).
- 35. Download the 'Current Results' csv file from the QIAcuity Software Suite.
- 10.2.1 Lambda Normalization
 - 36. The antibody concentration in each well of a QIAcuity Nanoplate 26k 24-well is reflected by the lambda value. The lambda values (in particular the distribution of the lambda values of one sample group) serves as a marker for pipetting precision. Normalizing the raw couplex counts based on their lambda values can compensate for pipetting inaccuracy.

Please note that lambda normalization can not completely compensate for pipetting inaccuracy and not for all cases of pipetting inaccuracy, therefore precise pipetting is important to obtain reliable results. 37. Calculate the lambda (λ) (average number of detected targets per partition) of each replicate (from each well of a QIAcuity Nanoplate 26k 24-well) using the 'Current Results' file.

Note: The lambda normalization for the raw couplexes can only be applied for one color. By default, we advise to apply the lambda normalization of the first color of the couplex pair as given by the AMULATOR export (column A) as 'type' informs about the color combination of the couplex. For example 'GY' stands for Green-Yellow. In this case, the lambda normalization will be performed using the lambda values of the green channel.

 $\lambda = ln(\frac{valid partitions}{negative partitions})$

Example: Calculate the lambda values for S#2 (wells A2-D2), S#3 (wells E2-H2), S#4 (A3-D3) and S#5 (wells E3-H3). As an example, the lambda normalization and ABC correction for well A2 is performed. The number of valid partitions for well A2 is 25,427. The number of negative partitions for the green channel is 21,025. The lambda value for this sample for the green channel is 0.190.

38. The Lambda Normalization Factor (NF_{λ}) is calculated for each sample (always for the color given first in the AMULATOR).

$$NF_{\lambda} = \frac{\lambda_{replicate} - avg(\lambda_{sample})}{avg(\lambda_{sample})}$$

 NF_{λ} : Lambda Normalization Factor for one replicate in the sample $\lambda_{replicate}$: Lambda value of the replicate

 $avg(\lambda_{sample})$: Average of lambda values of the sample

Example: The lambda value for the replicate in well A2 for the green channel is 0.190. The lambda values of the three other replicates for S#3 of wells B2, C2 and D2 for the green channel are 0.168, 0.191 and 0.156. The average of the lambda values is 0.176. The Lambda Normalization Factor for the sample in well E2 for the green channel is 0.078.

39. To calculate the lambda normalized couplex counts, first, the average raw couplex count of a sample has to be calculated ($avg(\xi_{sample})$). Then, the absolute value of the average has to be calculated ($|avg(\xi_{sample})|$).

Example: The raw couplex count for this replicate is 489. The raw couplex counts for the replicates in well B2, C2 and D2 are 470, 526 and 498. The $|avg(\xi_{sample})|$ is 495.75.

40. Calculate the lambda normalized couplex counts (ξ_{λ}) for each replicate using the raw couplex counts from the AMULATOR export.

Please note that the calculation of the lambda normalized couplexes does not influence the mean of the couplex count of a sample group but reduces the standard deviation.

$$\xi_{\lambda} = \xi - (\left| avg(\xi_{sample}) \right| \times NF_{\lambda})$$

 $\boldsymbol{\xi}_{\boldsymbol{\lambda}}$: Lambda normalized couplex count of one replicate

 ξ : Raw couplex count of the replicate

 $|avg(\xi_{sample})|$: Absolute value of the average raw couplex counts of the sample

 NF_{λ} : Lambda Normalization Factor for the replicate

Example: The Lambda Normalization Factor for the replicate in well A2 for the green channel is 0.078. The raw couplex count for this sample is 489. $|avg(\xi_{sample})|$ is 495.75. The lambda normalized couplex count for well A2 is 450.183.

10.2.2 ABC Correction

41. In the next step the lambda normalized couplex counts will be corrected to the ABC. In an ABC reaction, containing only the corresponding antibody mix without the sample, the antibodies colocalize in the compartments only according to Poisson statistics. Therefore, the couplex count in the ABC should be zero. However, due to offsets in the dPCR (e.g. signal dropouts or incorrect clustering) the ABC value may deviate from zero.

Note: ABC correction should only be applied when the ξ_{ABC} values derive statistically significant from zero, which is checked with a t-test.

Example: The raw couplex counts for the ABC control sample are -50, -45 and -85. The values derive statistically significant from zero (p-value: 0.0413).

42. First, the ABC Correction Factor (CF_A) for an experiment is calculated as the average of the raw couplex count of the ABC control sample.

 $CF_{A} = avg(\xi_{ABC})$

 ξ_{ABC} : Raw couplex counts of ABC sample

CF_A: ABC Correction Factor

Example: The raw couplex counts for the ABC control sample are -50, -45 and -85. The ABC Correction Factor, the average value of the ABC couplex counts, is -60.

43. The ABC corrected couplex counts (ξ_A) for each replicate is calculated using the lambda normalized couplex counts (see *10.2.1 step 41*) and the ABC Correction Factor.

$$\xi_A = \xi_\lambda - CF_A$$

- $\boldsymbol{\xi}_{\!A}\!\!:$ ABC corrected and lambda normalized couplex count of one replicate
- $\boldsymbol{\xi}_{_{\!\boldsymbol{\lambda}}}$: Lambda normalized couplex count of the replicate

*CF*₄: Average couplex count of ABC control sample replicates

Example: The lambda normalized couplex count for the replicate in well A2 is 450.183. The ABC corrected, lambda normalized couplex count (ξ_A) for the replicate in well A2 is 510.183.

10.3 Relative Quantification (RQ)

to S#2 and S#5.

- 44. For relative quantification ensure to use couplex counts from the low antigen side of the curve within the linear range. RQ is usually applied to relatively comparing target amounts in different sample types. In the case of RQ demonstration in this PICO Protein Detection Trial Kit, RQ is applied on samples within one dilution series which can also be used in other experimental setups to verify the data quality of a dilution series. At the low antigen side of the CLC curve, the raw couplex count is decreasing for increasing dilution of the sample and therefore the measured target (see 06 Scientific Background of a CLC Experiment). In the
- 45. First, test if the lambda-normalized and ABC-corrected couplex counts (ξ_A) are within the measurable range of the assay and not below the detection limit by testing if they ξ_A values of a sample are statistically significant different from zero.

experimental setup of the PICO Protein Detection Trial Kit, this corresponds

Example: The ξ_A values for S#2 (510.18, 552.15, 545.35, 615.60), S#3 (382.45, 309.67, 268.40, 278.14), S#4 (176.25, 205.98, 117.74, 220.02) and S#5 (-30.40, 21.56, -1.63, 94.46) are used to test if the measured dilutions of the recombinant human ErbB2/HER2 Protein are above the detection limit. The measured data of the samples S#2, S#3 and S#4 are statistically significant from zero, as tested with a two-sample t-test (p-values: S#2: 0.0001, S#3: 0.0012, S#4: 0.0046). In the case of S#5 (240,000-fold diluted sample), the sample input is below the detection limit (p-value: 0.1718) and will be excluded from the following RQ analysis.

46. In the case of the PICO Protein Detection Trial Kit, the applied dilution series of the sample and the known concentration of the recombinant human ErbB2/HER2 protein stock concentration (0.529 μ mol/l) offers the opportunity to evaluate the sensitivity of the PICO assay in this experimental setup. The values of the 24,000-fold diluted sample (S#4) are within the measurable range of the assay. Based on the setup of the binding reaction (2 μ l of sample is combined with 2 μ l of the ABX), the concentration of HER2 protein in the binding reaction is 48,000-fold diluted from the stock concentration, resulting in a concentration of 1.10 × 10⁻¹¹ M (equal to 11 pM).

Note that the theoretical limit of detection (based on analytical PICO models) is 4 couplexes and the theoretical limit of quantification is 10 couplexes. Due to the standard deviation, the experimental LOD is higher as demonstrated in this exemplary experiment. Based on this example, the average raw couplex count of S#4, with the lowest measurable protein concentration, the average couplex count is 12. Thus, one-twelfth of the couplex count measured in S#4 could theoretically be measured and quantified leading to a theoretical LOQ of 0.9 pM and one thirtieth of the couplex count could theoretically be measured, leading to a theoretical LOD of 0.37 pM.

47. The lambda-normalized and ABC-corrected couplex counts of the replicates (ξ_A) from the samples within the detectable range are normalized to the amount of recombinant human ErbB2/HER2 protein in the binding reaction (ξ_T). This is necessary to compare samples with different amounts of input.

Note that depending on the sample type, the input in the binding reaction can be a measure of target (in this case the amount or concentration of a recombinant protein) or a cell count in the binding reaction.

$$\xi_T = \frac{\xi_A}{T}$$

 ξ_r : ABC corrected and lambda normalized couplex count

of one replicate normalized to input of the binding reaction

 ξ_A : ABC corrected and lambda normalized couplex count of the replicate

T: Amount of input to the binding reaction

Example: The amount of recombinant human ErbB2/HER2 protein in the binding reaction of S#2, S#3 and S#4 is calculated based on the dilution series. The concentration of undiluted protein is 0.529 μ mol/l. According to *09.3 step 5*, each binding reaction contains 2 μ l of protein dilution. For S#2 a 6,000-fold dilution, for S#3 a 12,000-fold dilution and for S#4 a 24,000-fold dilution of the protein stock was performed. Thus, each binding reaction sample of S#2 contains 88.2 pmol, S#3 44.1 pmol and S#4 22 pmol. Exemplarily the ξ_A of the sample in well A2 belonging to S#2 is 510.18 resulting in a ξ_T of 5,785.48.

48. To evaluate if two data sets which are compared have statistically significant different expression of the target, the ξ_T values of both samples are tested for statistically significant differences. Choose an appropriate statistical test depending on your experimental setup. Based on the theoretical background of PICO, a normal distribution of the data can be assumed.

Example: Due to the experimental setup of the PICO Protein Detection Trial Kit, ξ_T values of the samples S#2, S#3 and S#4, all differently diluted samples from the same recombinant ErbB2/HER2 protein stock are compared. Thus, a paired t-test can be used. Further, we do not observe statistically significant differences between all samples (p-values: S#2/S#3: 0.424, S#2/S#4: 0.137, S#3/S#4: 0.378) as expected, because the data was already compensated for the different amount of input. (The ξ_T values for all samples can be calculated based on the ξ_A values given in the example of step 21 and the sample input given in example of step 23.)

49. Next, the fold change between to samples is calculated using the average ξ_T values of two samples.

 $FC = \frac{avg(\xi_T(sample 1))}{avg(\xi_T(group 2))}$

FC : Fold change of the ξ_r between both samples

 $avg(\xi_r(sample 1/sample 2): Average lambda normalized, ABC$

corrected and input normalized couplex counts of sample 1 or 2

Example: The fold change using the average ξ_T values of S#2 (5,836.11), S#3 (6,503.00) and S#4 (7,560.00) samples were calculated resulting in FC_{S#2/S#3} of 0.897, FC_{S#2/S#4} of 0.772, and FC_{S#3/S#4} of 0.860.

50. Transform the fold change values into log2 fold change values (results in $log_2(FC)$). This value indicates the relative difference of the analyzed target present in the different samples (meaning the differences in the initial undiluted sample, not the diluted sample used for RQ).

Example: The fold change values are transformed into log2 fold change values $(\log_2(FC)_{S\#2/S\#3}: -0.156, \log_2(FC)_{S\#2/S\#4}: -0.373, \log_2(FC)_{S\#3/S\#4}: -0.217)$. As expected, the values are close to zero, indicating no different expression between the samples.

51. The p-value can be transformed using $-\log_{10}(\text{p-value})$). The log2 fold change can be plotted against the transformed p-value in a volcano plot.

When RQ is performed to compare the amount of target in two samples, we recommend to average the log2FC values and the p-values.

Example: The p-values ((p-values: S#2/S#3: 0.4241, S#2/S#4: 0.1369, S#3/S#4: 0.3778)) are transformed into -log10 values (S#2/S#3: 0.3725, S#2/S#4: 0.8635, S#3/S#4: 0.4227). The average log2 fold change is -0.249 and the average -log10 p-values is 0.553. Thus, in case of the experimental setup of the PICO Protein Detection Trial Kit, the relative differences between S#2, S#3 and S#4, all differently diluted samples from the same recombinant ErbB2/HER2 protein stock, result, as expected, in no significant difference in protein amount.

11 Troubleshooting Guide

Consult the troubleshooting guide below to solve any problems that may arise. For more information, see the <u>Frequently Asked Questions</u> (FAQ). It is available at: <u>www.actome.de/resources/downloads</u>

Reach out to the Actome Customer Service (<u>info@actome.de</u>) if you have any questions about the protocol.

Troubleshooting				
Issue	Comments and Suggestions			
No couplexes or low numbers of couplexes detected				
Wrong default threshold of fluorescence intensity (RFU) was set in the QIAcuity software suite	Select '1D Scatter Plot' in analysis mode of QIAcuity software suite and adapt the thresholds.			
Error from AMULATOR				
Single wells of the nanoplates are empty	Delete the corresponding wells in the sample definition file as well as the according table line to create a continuous table.			

12 Ordering Information

PICO kits can be purchased directly from <u>shop.actome.de</u> or a quote can be requested from <u>sales@actome.de</u>. Supporting materials are available on <u>www.actome.de/resources/downloads</u> or can be requested from Actome's Customer Support (<u>info@actome.de</u>).

Ordering				
Product	Description	Cat.#		
This Product				
PICO Protein Detection Trial Kit	Trial kit for the detection of recombinant HER2 using trastuzumab and pertuzumab	PICO-000090		
Related PICO Products				
PICO Amplification Core Kit	dPCR detection for PICO assays (5 x 24 rxns)	PICO-000010		
PICO Probes	BL (HEX), P8 (FAM), N6 (TAMRA), O7 (ROX) probes for detection in dPCR (5 x 24 rxns)	PICO-000070 - 73		

